

Autoclave Maintenance




HMC 206FA in ETH Biolab (D92.2)

Daily Routine (after use)



1. wipe the chamber with a cloth
2. check the water level of the tank, refill when low. Use only RO / distilled water to refill
3. check that the rubber lid of the water tank is in place and allows air to pass through the airholes

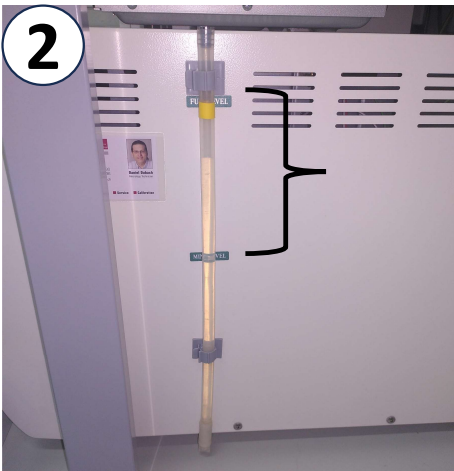


Weekly check-up


1. Take out the metal tray and clean it



2. check all openings in the water tank and clean/unclog them if necessary
3. Pull on the safety pin 3x to check function

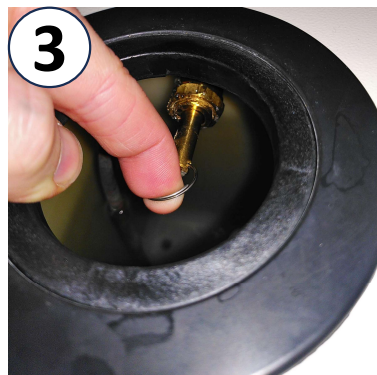

This is how the water tank should look like

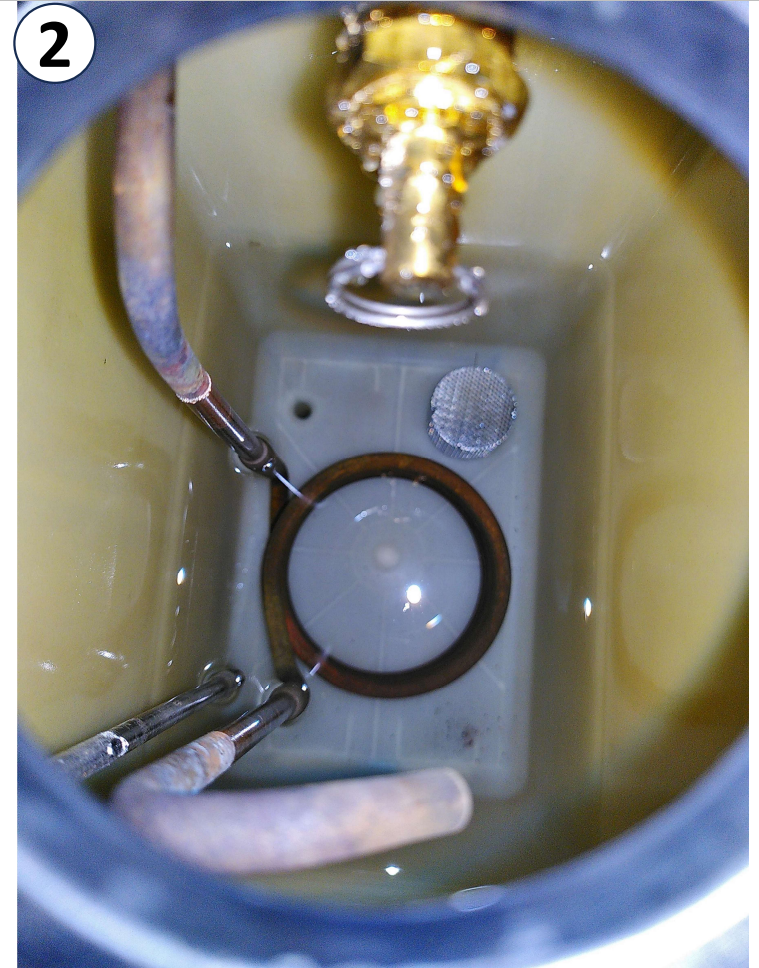

Monthly cleaning (1/3) clean the water tank as follows:



1. release the water tube
2. remove the metal cap for faster flow
3. lower the water tube and collect the water in a tray until tank is empty



Monthly cleaning (2/3)



Rinse the water tank by pouring water into the top

- 3x with 1.7 L cold water
- 3x with boiling hot water 1.7 L
- reinstall the water tube and refill the tank with RO water

Monthly cleaning (3/3)



1. Locate the filter at the back of the autoclave
2. Place a towel underneath to prevent spills
3. Unscrew the filter
4. Rinse it with water and clean out debris and screw the filter back in

